

Bel-Art Products
6 Industrial Road
Pequannock, NJ 07440
Tel: 973-694-0500 Fax: 800-545-9796

Refilling Eye Wash Bottles

1. Water only – change once a week
2. Water and Preservative – vary from preservative manufacturers, change every two months
3. Water and Eyesaline Concentrate change every four months
4. Pre mixed Eye Wash Solution every 6 months
5. Sterile Eye Wash Solution unopened until expiration date on bottle.

The above recommendations for solution replacement are based on information gathered from sources outside Bel-Art Products to help assist our customers. Bel-Art Products does not guarantee the accuracy of the information gathered.

Bel-Art Products Sterile Eye Wash solution (24879-0032) is not meant to be used as a replacement in refilling eyewash bottles, as the integrity of the Sterility of the solution is jeopardized once the bottle is opened. It is meant for use with our 24878-1032 and 24878-2032 Eye Wash Stations. Refer to expiration date on bottle for shelf life normally 3 years.

If using the Bel-Art Products Sterile Solution to refill bottles with eye cups, the solution can no longer be considered sterile once transferred into the those bottles. Replacement of solution once transferred should be 4 to 6 months.